



## IHR 320 Blaze Value Change

This procedure explains how to change the blaze values stored for a Horiba iHR320 Spectrometer. This assumes that the software has already been installed and that the specific spectrometer has been set up for use. If you have multiple units, you will need to select the particular unit in the Spectrometer drop down menu shown in step 1.


1. Open USBSpectrometer Control

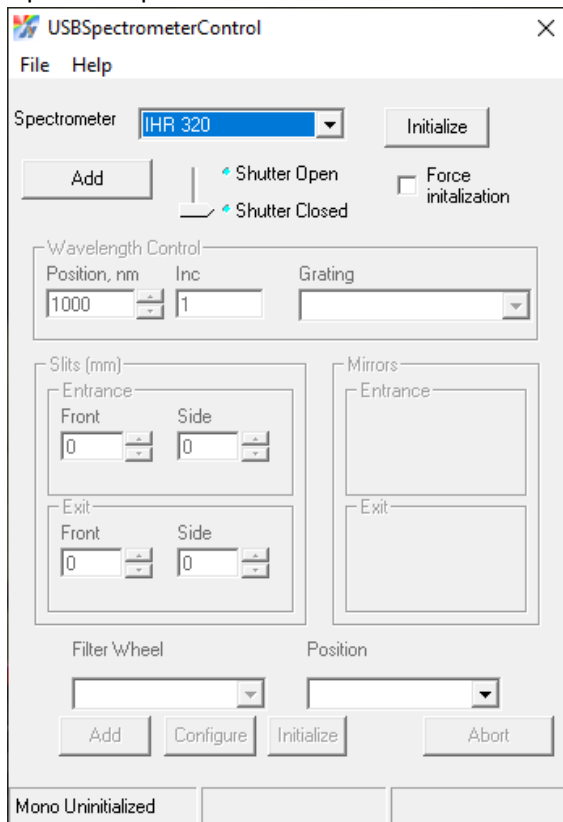


2. Select File Menu, then Configure

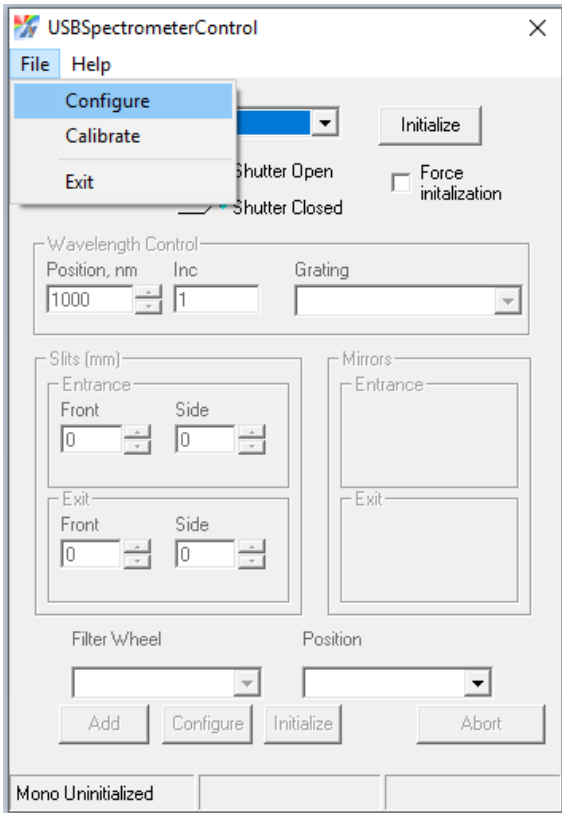


3. Select AccessoryConfig

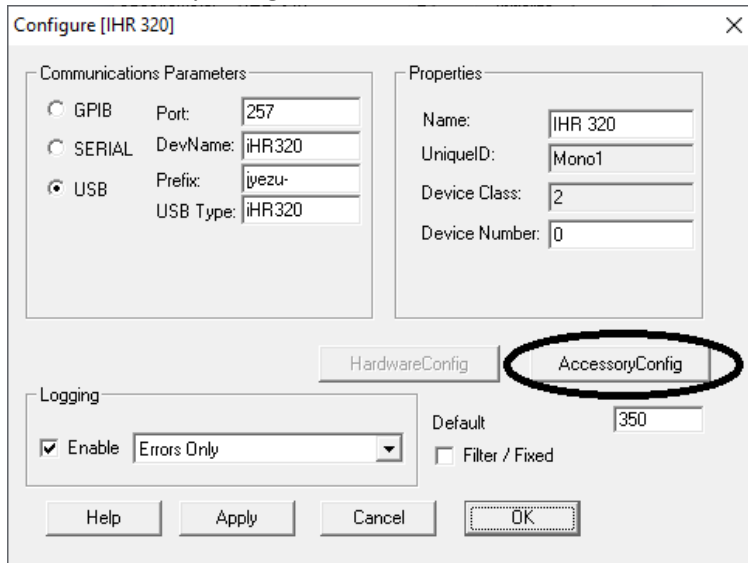

4. Change Blaze Number/Text values and click OK

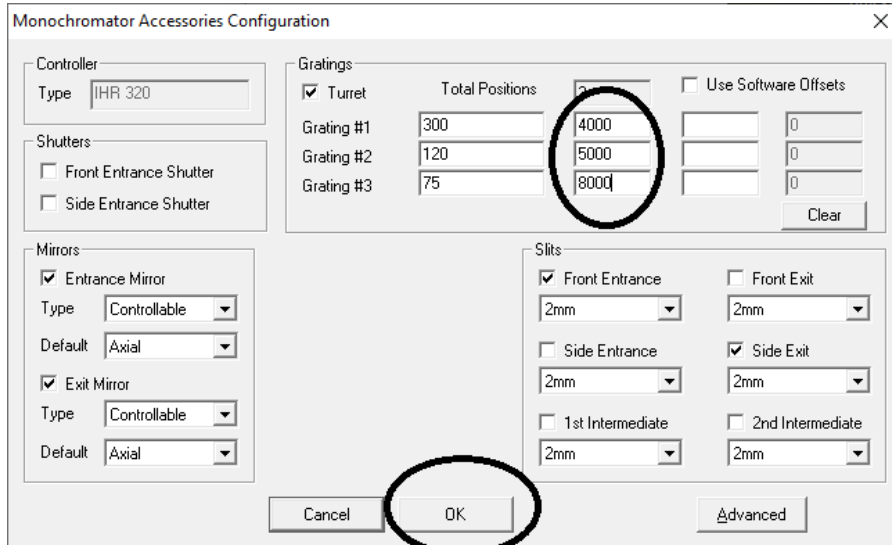

5. Click Apply, then OK.

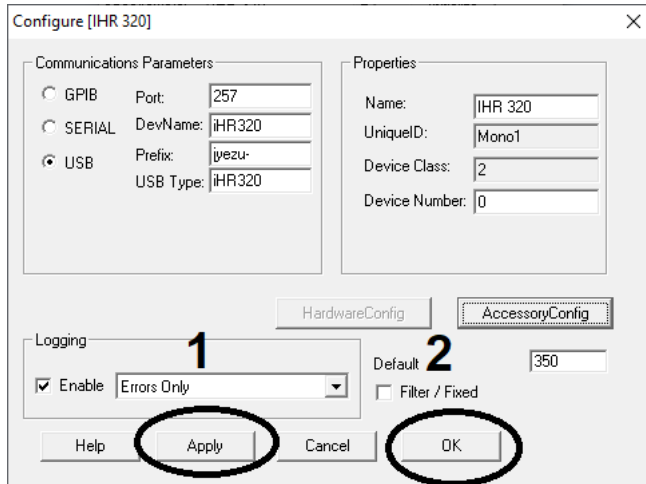

6. Reinitialize (Force Initialization) to use new values. Procedure is now complete.